



adolphe merkle institute
excellence in pure and applied nanoscience

UNIVERSITY
OF FRIBOURG
SWITZERLAND

Focused ion beam

Introduction

Version 1 – February 2022

FIB-SEM EDX



Quick summary

Start

- Setup the FIB-SEM: Load sample, pump chamber, eucentric height, navCam, ...
- Get an image in the SEM
- Stop the chamber view camera
- Assure both PCs have the user logged in ('user')

Counts

- Insert the detector and start the Pathfinder software
- Assure the SEM is acquiring
- Spectrum > start spectrum dropdown > high resolution > advanced arrow
- setup the SEM to have stores > 2000 and a dead time < 40-50% (ideal: 20-35%)

Spectrum

- Collect a spectrum
- Identify the elements

Data

- PointID
- Spectral imaging
- Linescan

End

- Export the data from the database
- retract the detector
- Extract your samples, set the FIB-SEM to standby
- Cleanup the desk

Universal rules

Rule 1: don't touch a control if you are not sure of the outcome of that action

Rule 2: never, ever force anything beyond finger strength

Rule 3: wear gloves when touching anything that goes into the chamber

Rule 4: if in doubt, ask for help

Demonstration: Ending your session

Prerequisites:

Data recorded and exported

Ready to extract your data from the database

Experiment: end the EDX session

1. Switch on the chamber view
2. Retract the detector entirely (must not be visible anymore in chamber view)
3. Close the pathfinder software
4. Stop the SEM beam and put the FIB-SEM in standby mode

Demonstration: setup the microscope6

Demonstration: setup the beam7

Demonstration: Start the pathfinder software.....8

Demonstration: optimizing X-ray counts.....10

Demonstration: A word on the time constant.....12

Demonstration: Record a spectrum13

Demonstration: Identify using the periodic system14

Demonstration: Point ID mode.....16

Demonstration: Spectral imaging.....17

Demonstration: Adjusting the spectral maps.....18

Demonstration: Extracting spectra from maps20

Demonstration: Map quantification22

Demonstration: Multivariate statistical analysis with COMPASS.....23

Demonstration: Phase maps.....24

Demonstration: X-ray topography.....25

Demonstration: Line scan26

Demonstration: Saving your data27

Demonstration: Ending your session28

Demonstration: setup the microscope

Prerequisites:
Functional FIB-SEM

Starting your session

You do not need the FIB beam; the Ga ion beam is destructive and does not produce X-rays.


Experiment

Make sure you are logged in with username 'user' on the FIB-SEM microscope PC

Setup the basics at the FIB-SEM/SEM imaging

- Load your sample
- Home the stage
- Put the sample in eucentric height / 7 mm
- Make a navcam picture
- Use the ETD for imaging

Assure that no large sample is blocking the way to right ride of the chamber (in chamber view). Below: suboptimal: the sample on the right might block the path

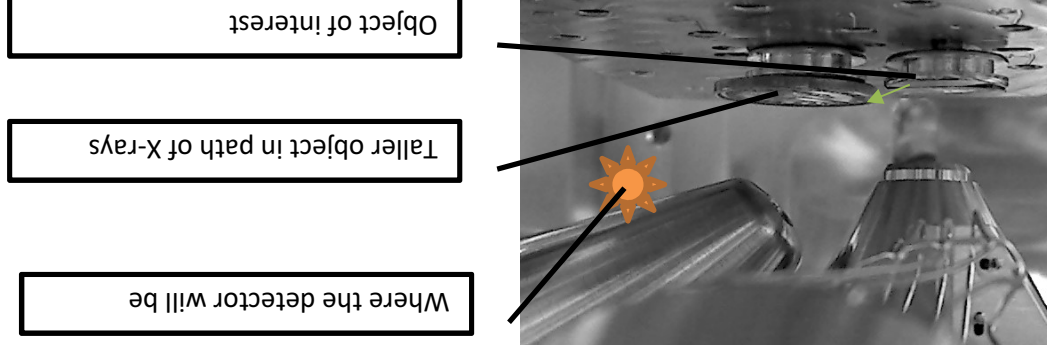

Demonstration: Saving your data

Prerequisites:
Data recorded

Ready to extract your data from the database

The data is automatically saved in the database. To extract the data from the database:

1. Export full field spectrum

- Look into your folder where you saved your database.
- There will be an **.emsa file** for every full spectrum you recorded⁵.

2. Point ID spectra

- Look into your folder where you saved your database.
- There will be a **folder** with the base name of your recording.
- Inside the folder you will find a **.psmsa file**. This is equivalent to the **.emsa files**.

3. Extract spectral imaging maps

- Save the maps by file > save map files.
- You will be asked to give the map a type (e.g. counts)
- A new entry will appear in the data list (e.g. filename.counts).
- The maps of all elements can be found in the database root folder as tif files (using the color code as setup in Pathfinder).

3. Extract line scans

- Select your line scan
- Save as CSV or tiff by file > import / export

⁵.emsa files are text files with the spectrum and metadata.

Demonstration: Line scan

- Prerequisites:
 Detector inserted
 Stores / s optimized

Record a line scan at high signal to noise

The functionality of the linescan is similar to the line scan extraction by spectral imaging. However, by only scanning the points along the line (and not the entire map), a longer integration time can be achieved and thus a much better signal to noise ratio.

Experiment: carry out a line scan

- In the microanalysis tab: click Linescan
- Similar to the spectral imaging, record a SEM Image and then record a linescan.
- Adjust the resolution of the scan if needed



Bug alert: You must save your file (File > close project) and reopen the database in order to adjust something to the output.

Demonstration: setup the beam

- Prerequisites:
 Sample(s) loaded
 SEM bean on


Starting your session

Over-voltage

The choice of the high voltage must be seen in the light of the elements you want to detect. As a rule of thumb:

$$HV \geq \frac{E_{x-ray}}{0.5 \text{ to } 0.7}$$

Or at 30 kV (the maximum on the FIB-SEM), you can detect up to 21 keV edges. Use the Slider tool to extract the information. For example, it Au is the goal:

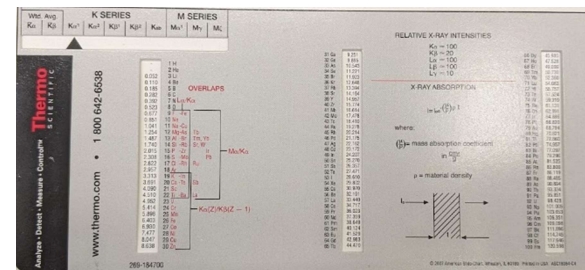

- The K_{α} edge of Au is at 68.794 keV (= undetectable)
- The M_{α} edge of Au is at 2.123 keV (= detectable)
- The L_{α} edge of Au is at 9.711 keV (= detectable if HV > 13.9 kV).

Therefore, 15-20 kV is a good setting to detect Au.

Beam current

The beam current can be set to standard values, e.g 0.4-0.8 nA.

Magnification

Do not zoom in too much to start. 500-1000X

Demonstration: X-ray topography

Prerequisites:

Spectral maps recorded

3D graphs of the X-ray distribution


Experiment: X-ray topography plots



X-Ray Topography



- Select surface in the cursor tab
- The workspace changes. A large white field appears on the top left.
- Select an element in the maps
- Choose your settings in the X-ray topography tab

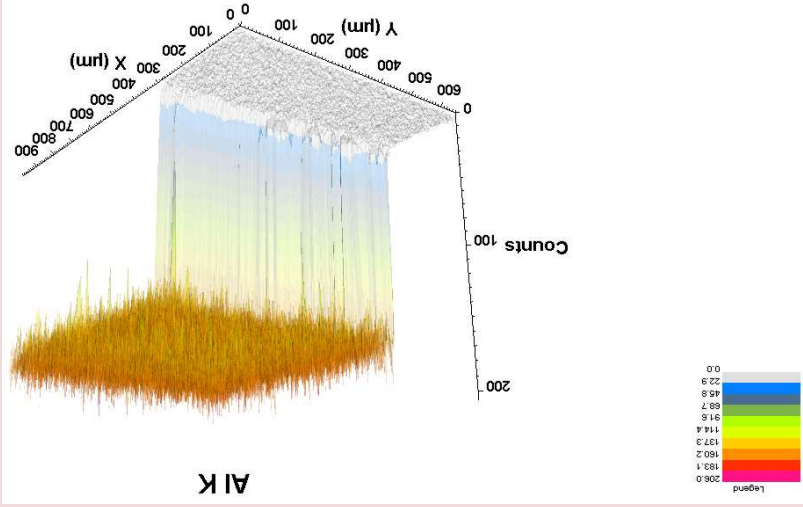

Demonstration: Start the pathfinder software

Prerequisites:

Proper image under the SEM

Logged in under the username 'user' on the Suppoer (EDX) PC

It is absolutely crucial that both PC are logged in with the same username: 'user'

Before you start: On the right side of the table, check the LED's on the SSD controller. All four LED's must be green.



Experiment: Insert the detector

- Insert the Detector using the EDX switch on the table. Press insert for 2-3 seconds until you see the detector extended into the chamber view. The LED on the switch will turn red.

EDX detector retracted



EDX detector inserted



Demonstration: Phase maps

Prerequisites:
Spectral maps recorded

Extract compositionally distinct binary phases

Experiment: use Phases

In Map processing, find phases. You get a list of the elements (uncalculated phases)

Phase map type

Maximum intensity:	Uses MI in each pixel to create the phase maps
X Phase:	proprietary algorithm
Auto	No user input needed
Manual	Allows to set threshold

Experiment: start the Pathfinder software and login onto a database


- Start the Pathfinder (icon on the desktop) on the EDX PC



- A splash screen is visible for 10-15 seconds.

- You are asked to use a database. Use an existing one or create your own in your user folder.

- **You do not have to actively start a cooling**

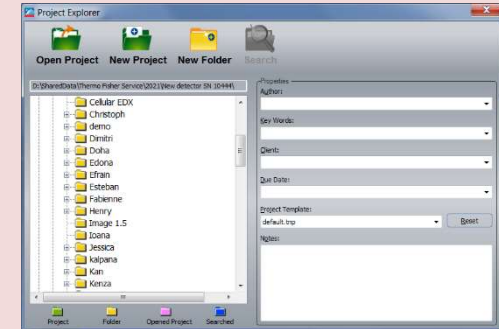

Demonstration: optimizing X-ray counts

Prerequisites:


Detector inserted

This important setting is very hidden in the software

Before you start recording data, you should optimize the X-ray counts.


Experiment: finding the X-ray counts

- On the top right, in the microanalysis tab, choose spectrum:

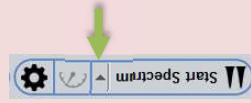

- The click the down arrow next to 'Start spectrum' (Do not start a spectrum!) and choose any setting (e.g. High resolution).

- Then, in the top right of the experiment setup, click the green down arrow

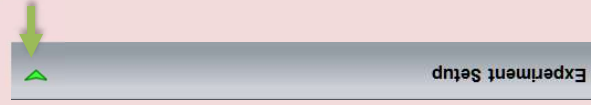

- Now, the counts and stores are at the bottom of the Experiment setup.



Experiment: optimizing the X-ray counts

1. Switch off the chamber view, as this also produces X-rays. Click the lower right quadrant and press F6. Make sure the topleft corner has this symbol:

Demonstration: Multivariate statistical analysis⁴ with COMPASS

Prerequisites:

Spectral maps recorded

Extract compositionally distinct components without user input

Experiment: use COMPASS

Method

Area: weight by nearest neighbor, Recommended for most sample types

Spectral: No NN. Intended for samples with many small phases

Background


Using data: for data with sufficiently high signal to noise

Using model: for noisy data

Number of components

0 (auto) recommended as pilot

1-15 optimization options

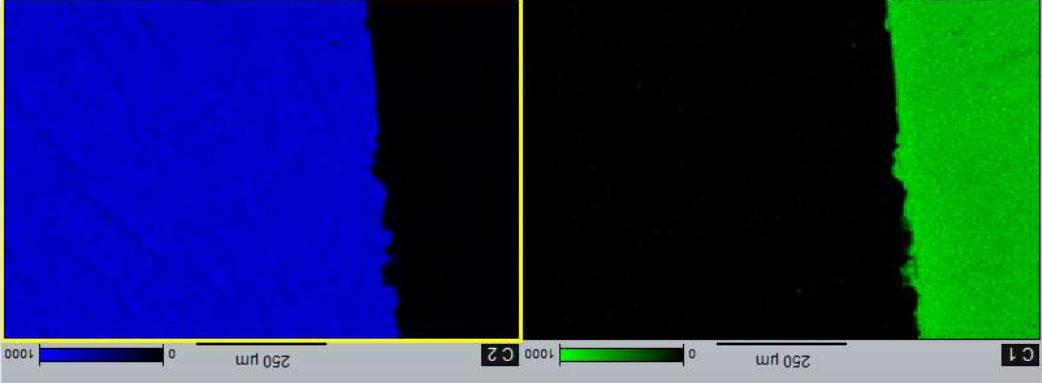

⁴ Based on principal component analysis


Demonstration: Map quantification

Prerequisites:
Spectral maps recorded

Converting counts into atomic %

Experiment: convert counts to atomic %

- Select QUANT in map processing
- Deselect 'threshold' to assure you process all maps
- Set the endpoints you want to achieve:
Atomic %, weight %, ...
- time you want to invest (with 256px maps, it is a matter of seconds)
- Kernel will apply a mean filter, increasing the signal to noise at the cost of resolution.
- Click process

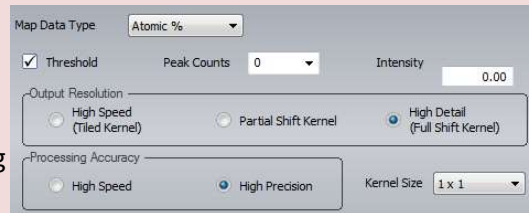


The result are maps with a dynamic range of 100 (from 0% to 100% atomic or weight %)

Intensity cursor




x: 147 y: 67
 Grey: 1.34%
 C K: 0.00%
 O K: 2.24%
 F K: 0.00%
 Mg K: 5.37%
 Al K: 92.08%
 Cl K: 0.00%
 Co L: 0.00%
 Cu L: 0.00%
 Pd L: 0.31%
 Au M: 0.00%

Now, use intensity cursor ('+ cursor') to output the atomic % for each element at a specific map (xy) position.

2. Start the SEM scanning (i.e. there is NOT the symbol  on the top right quadrant)¹
 - Image size: e.g. 768 x 512
 - Scan speed for good imaging: 5-10 μs/pixel (depends on the image size)
 - Detector: ETD
3. Adjust the Beam current until you have
 - high stores count (> 2000, better > 10 000)
 - not too high dead times (15 > dead time < 40)

Dead time

- High dead times means inefficient collection
 - High dead times cause artefacts, such as higher sum peaks
- A high dead time means the detector is swamped with X-rays while the collection

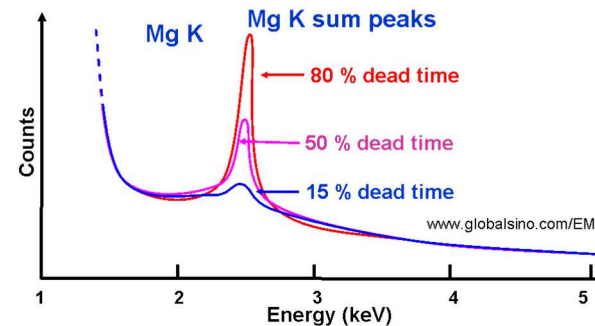

- However, lower dead times may mean lower stores / s

Example of a good setting

Detects /s:	18800	<div style="width: 100%; height: 10px; background-color: green;"></div>
Stores /s:	13592	<div style="width: 100%; height: 10px; background-color: green;"></div>
Dead Time %:	23	<div style="width: 100%; height: 10px; background-color: green;"></div>

¹ At some point, the software can activate the scanheads, but not at the first scan.

Demonstration: A word on the time constant

Prerequisites:

Detector inserted

Stores / s optimized

The time constant affects the quality of your detection

X-rays generated from a specimen are subjected to dispersion according to their energy. during acquisition, averaging of this noise is performed to obtain accurate values.

The time constant is a time index to define the noise collection. Changing the time constant, the averaging time for noise collection is changed.


High time constant → good average

high-energy resolution EDS spectrum
higher sum peaks, low throughput

Low time constant → poor average

lower dead time, efficient imaging,
high throughput
More noise = lower spectral resolution

Appears on the image.
- Adjust the line (grab the edges)
- On the right: linescan processing becomes available. Adjust if needed.

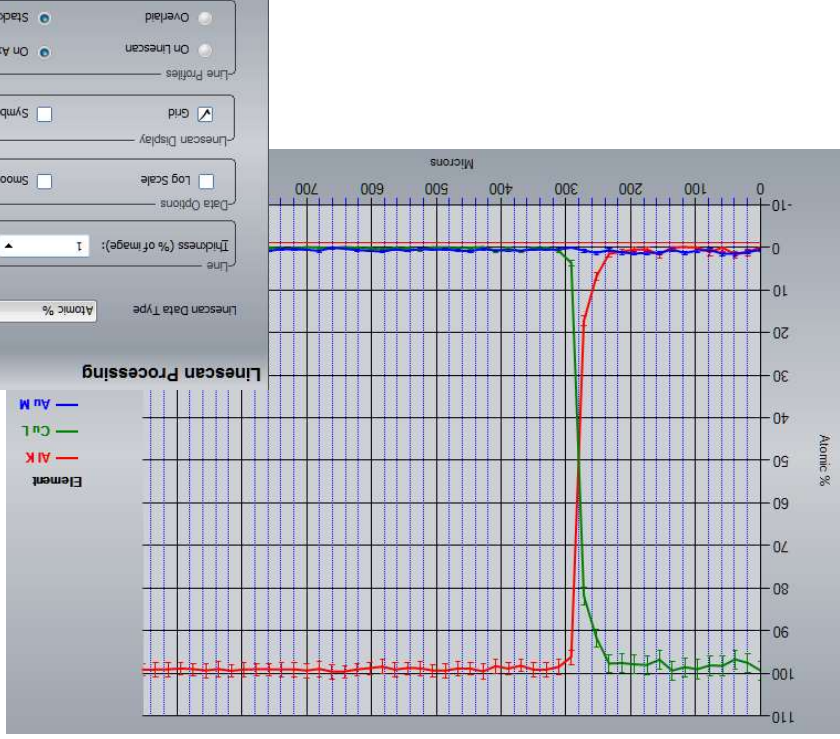


Demonstration: Extracting spectra from maps

Prerequisites:
Spectral maps recorded

Browse through the data in the spectral maps

1. Extract data from points or region of interests

In the bar under the SEM image, select a cursor type (rectangle, spot, ...) and select a region in the SEM/Overlay image. The extracted spectrum will show up in the spectrum pane (make sure intensity cursor '+ cursor' is switched off).

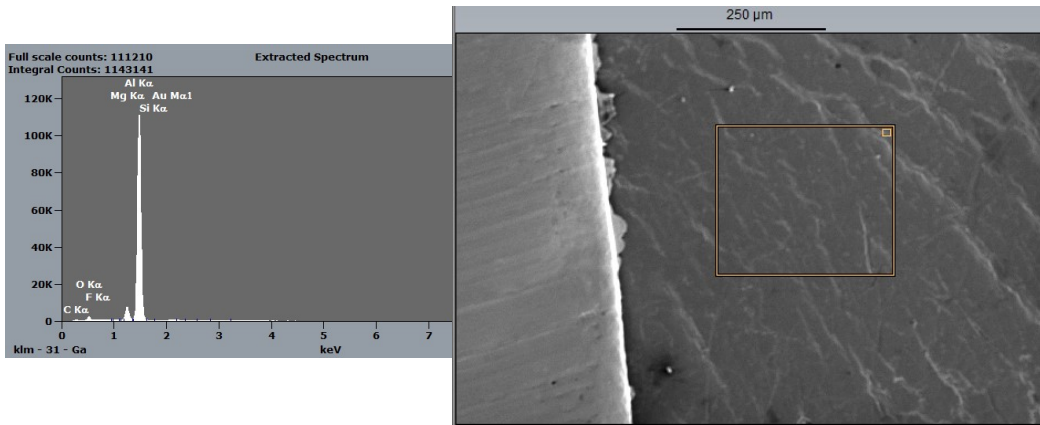


The spectral data is shown either as:

- Area integral Total X-ray counts from the selected area (default)
- Trace element search Maximum counts for each eV channel (useful for finding Small concentrations of an element)

2. Extract line plots

- Select 'Line'. An orange line overlay

D Vanhecke | Adolphe Merkle Institute

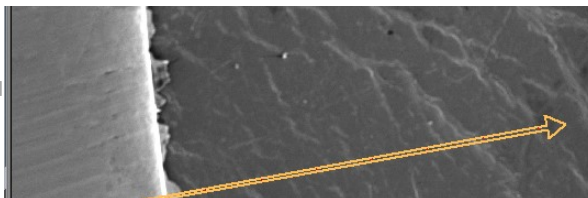

Demonstration: Record a spectrum

Prerequisites:
Detector inserted
Stores / s optimized

Record a spectrum of the entire field of view

The 'Experiment' control panel shows the following settings:

- Termination Criteria:** Live Time Limit (s): 30, Max Peak Counts: 10000
- Energy Range:** Low Energy Cutoff (eV): 100, High Energy Cutoff (keV): Auto
- Time Constant:** Rate 1, Shaping time 71400 cps (projected max throughput)

Annotations in the image:

- Stop the acquisition after e.g. 30 seconds, or a peak reaches e.g. 10 000 counts
- Cut off artefacts/noise below 100 eV. Leave HEC to Auto
- High throughput or high spectral resolution

Buttons: Click start spectrum, Start Spectrum, [Settings icon]

Result: Spectrum processing

- Assure that Auto ID is on & click the green arrow
- There are some options for the Auto ID
- Click process to repeat the analysis²



The 'Spectrum Processing' control panel shows the following settings:

- Ident Setup:** [Ident Setup icon]
- Quant Setup:** [Quant Setup icon]
- Auto ID:** [Auto ID checkbox]
- Peak Removal:** [Sum Peak Removal checkbox], [Escape Peak Removal checkbox]
- Ident Sensitivity:** 5
- Overtoltage:** 1.5

² Process will only start the identification of the peaks. Clicking Identify button in the bottom left will do the identification and the quantification.

Demonstration: Identify using the periodic system



Prerequisites:
Spectrum recorded

Record a spectrum of the entire field of view

KLM lines


With no auto ID run (all peaks unknown), left click on an element and the peaks will be overlaid on the spectrum.

Click left/right on search to go through all elements.

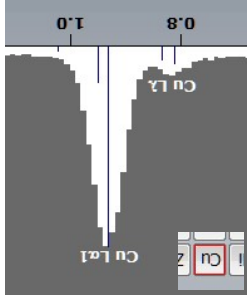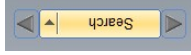

Select the peak element

If the element was detected, double-click to see other available elements.

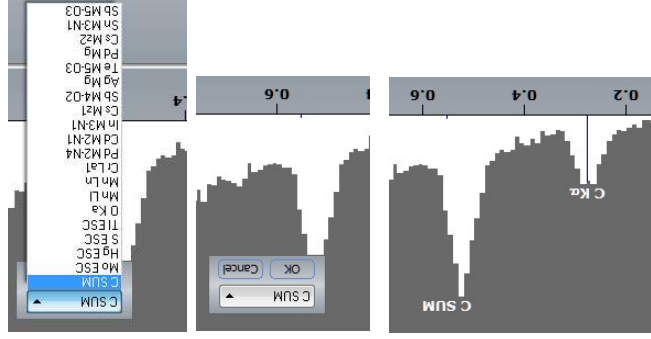


Compare with a synthetic curve

Comparing with a theoretical curve allows you to see if you missed any elements

- Filters will create severe artefacts in data with low dynamic range. Increase dynamic range by recording longer.
- In high dynamic range data, filters are not advisable.

The software allows to improve noise by filters (especially the smoothing filters).

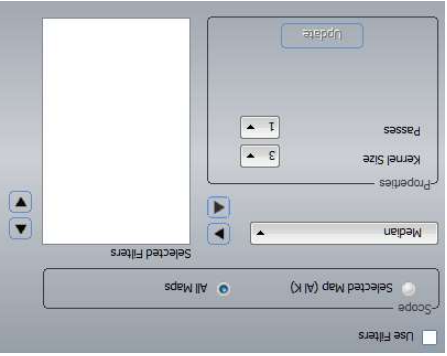

When you use filters, make sure you know what you are doing!

4. Change transparency of the overlays



- In Image settings, choose Electron

image.

- BUG ALERT: Click image filters and then again on image settings!!

- You can now change the

transparency of the overlays

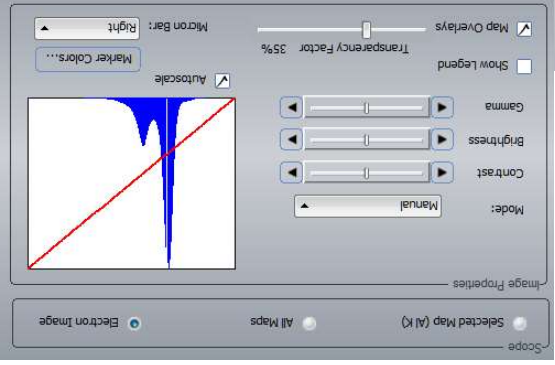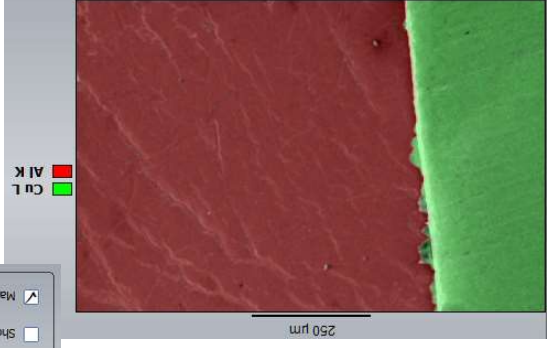

Demonstration: Adjusting the spectral maps

Prerequisites:
Spectral maps recorded


Change the visuals of the maps

1. Change the selected elements

Use the periodic table, right-click on the wrong element and choose 'excluded' to remove elements.

2. Change the false color

- Left click the map (a yellow border appears).
- Choose image settings
- Choose Selected Map (in this case Al K)
- **BUG ALERT: Click Image filters and then again on Image settings!!**
- Now you can change the colors


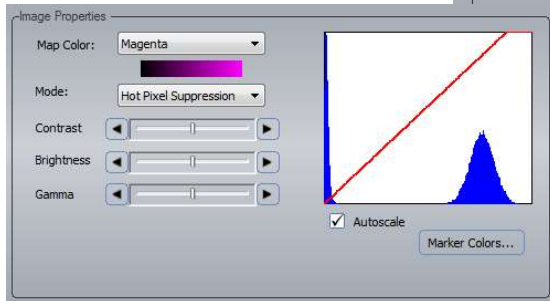


3. Image filters

Select your data from the bottom right and click the compare icon



- Leave normalization to none, but click the spectra check:
- 3 new entries become available in the list:
- click Synthetic spectrum
- the goodness of fit is shown by hovering over the quantification pane title. The lower the better.

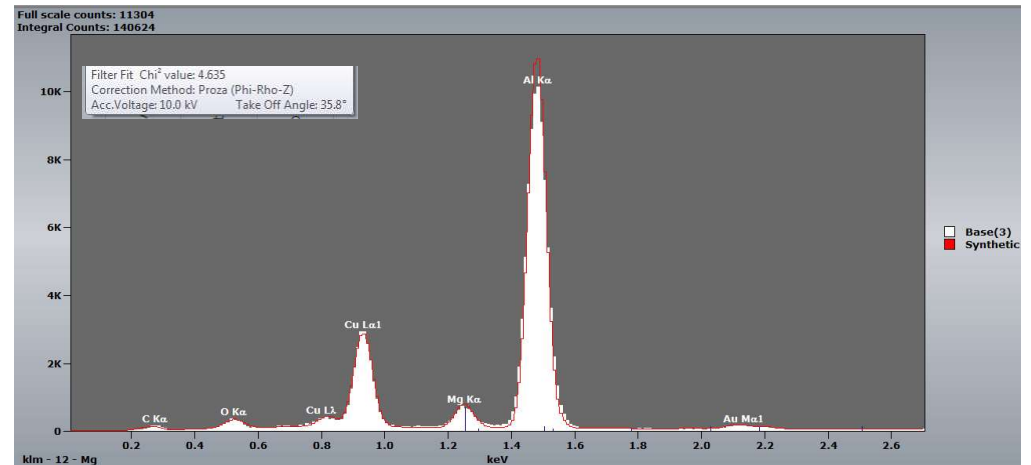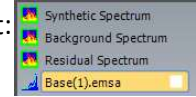

Demonstration: Spectral imaging

Prerequisites:

Detector inserted

Stores / s optimized

Record an elemental map

Experiment: acquire a map


1. Acquire an SEM image (see before)

2. Setup the variables. Map scans can take a very long time, and map size is a key setting affecting the time. Do not overdo the map size. Usually, 256px is a good setting.

- *Frame time*: Time scan 1 frame

- *Number of frames*: repetitions

3. Start the scan. If the autoid is on (standard), the maps of the detected elements will be shown in the central pane.

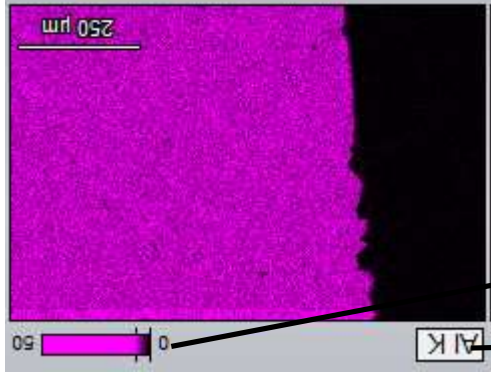

Element + overlay toggle

Min and max intensity

It is advisable to have at least 100 values in the scan (this is the dynamic range). More is better.

Demonstration: Point ID mode

Prerequisites:

Detector inserted

Stores / s optimized

Record a spectrum of the entire field of view

Experiment: get a point spectrum

1. Record an SEM image: Click Get Image after setting up the imaging details.

2. In the SEM image pane, select the geometry (point, ...)

And draw or place your

ROI. If instant is selected

The scan will start

Immediately. Otherwise,

Click start spectrum

