

Focused ion beam

Introduction

Version 10 – May 2025



ThermoFischer SCIOS 2 Introduction – Page 3

Universal rules

Rule 1: don't touch a control if you are not sure of the outcome of that action

Rule 2: never, ever force anything beyond finger strength

Rule 3: wear gloves when touching anything that goes into the chamber

Rule 4: if in doubt, ask for help

Agenda

Agenda	5
Demonstration: xT microscope server	6
Demonstration: Login onto the FIB software	7
Demonstration: System status	8
Demonstration: Loading a sample	9
Demonstration: Beam control	13
Demonstration: The inside of the e-beam	15
Demonstration: First glimpse of your sample	17
Demonstration: Eucentric height	19
Demonstration: Nav Cam photo	21
Demonstration: Basic SEM imaging	22
Demonstration: Beam and lens parameters	29
Demonstration: Imaging settings	30
Demonstration: Image presets	32
Demonstration: Align Feature / scan rotation	33
Demonstration: Column cases	34
Demonstration: Detectors	36
Demonstration: Direct adjustments – Source tilt	39
Demonstration: Beam deceleration	40
Demonstration: Finishing your session	42

Prerequisites: MPC (microscope PC) and SPC (support PC) switched on Login: User (password: user) Start the server Experiment Find the tiny view server bar at the top right of the central screen: 83 Start Stop Hide UI Stop UI X The xT server must run in order to use any FIB function. It runs when you arrive at the FIB. YOU NEVER LOG OUT, YOU NEVER SWITCH OFF THE PC's not enabled = server is running Start Stop enables = server is running Hide UI hides the user interface (will become start UI when switched off) Stop UI closes the user interface If the user interface is not active, start it by clicking "Start UI"

Notes:

Demonstration: Login onto the FIB software

Prerequisites:

Running xT server Running UI

Action:

Login using your FIB account

Load the personal settings and history of the user

Experiment

After startup of the User interface (UI), a username and password are requested.

	FEI microscope
- D	vanhecke
	password
	Log On

Username: your last name, with first letter capital, no accents, umlauts, etc. Password: your first name (first letter capital but no accents, umlauts, etc.)

Click logon

Demonstration: System status

Prerequisites:

Running xT server Running UI and successfully logged on

Notice the application status

A separate window will appear (usually on the rightmost screen): the application status, which will contain messages from the system. These will contain any errors that occurred.



You can close the application status window. You will see new messages in the bottom bar of the UI. Clicking the message button will reopen the application status window.

If the system behaves out of the ordinary, this is the first place to look for error messages.

Demonstration: Loading a sample

Prerequisites:

Running xT server Running UI and successfully logged on

Action:

Load samples

Vent the chamber. Understand the height prerequisites of the load samples

Click the sample exchange button



on the top left of the central screen.



Experiment: Venting the chamber

Under Chamber, click vent. A venting confirmation will ask you to confirm the 3 minutes venting cycle. Click vent and wait 3 minutes (the green bar at the bottom shows the progress).

The multipurpose holder



In most slots, the stubs can be mounted without the need for a screwdriver.

Experiment: mounting the samples

- Gently open the stage
- Place the stubs in an empty slot on the multipurpose holder.
- Make sure that you place:
 - at least 1 stub / sample in the central position (#9)
 - the tallest samples on the left,
 - the lowest on the right (this avoids pole touches while tilting).
- Gently close the chamber. There is no lock.

The Vacuum system

Holder: Always have the Multipurpose holder selected. Never change this setting.

Chamber:

- Sample-cleaning: starts a in-chamber plasma cleaning of the stage/samples. Can help with the conductivity of notoriously badly conducting materials

Never use the chamber plasma cleaning with the EDS detector inserted!

- The take Nav-Ca Photos does not make sense because the eucentric height has not been set yet.

Tab Vacuum

The value mentioned under TMP1 is critical for the pumping of the chamber (see below)



D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland

Experiment: pump the chamber

- In the sample exchange window. In summary:

- Sample cleaning (optional, not with an inserted EDS detector)
- Do not select 'Take Nav-cam Photo',
- Do not change the root folder settings.
- Make sure multipurpose holder is selected

- Switch to the vacuum tab and locate the TMP1 reading (bottom right, under 'Gauges'). It should have the value 0%.

Keep the door firmly closed by leaning against the chamber door with your arm. This is extremely important!¹

- Click the Pump button in the exchange window.

- Keep leaning against the door / holding it tight until you see the TMP1 value increasing to at least 5% of the TMP1 (you stop hearing the rotation pump in the back). Now you can take back your seat.

- While waiting for the vacuum, home the stage: *Stage* > home stage.

After pumping, you should reach a chamber pressure of about $1 \cdot 10^{-2}$ Pa. This is not yet sufficient for a focused ion beam ($1 \cdot 10^{-4}$ Pa is needed), but it allows e-beam imaging. Pumping will go on in the back and by the time you have setup the e-beam, the necessary vacuum for the ion beam will be achieved.

¹ If the door is not properly sealed, an overheat safety will halt the pump. The pump (and the system) can only be reset by a technician. Hence, not sealing the door will mean the end of your session and about 1 week of downtime for all other users.

Demonstration: Beam control

Prerequisites:

Sample loaded Pressure at $1 \cdot 10^{-2}$ Pa or lower

Starting the e-beam

On the right side of the central screen, an icon-based menu is available:

Beam control	Vacuum		1
Navigation	Pump Sample Cleaning	Vent	
Detectors	System		
Patterning	Wake Up Column	Sleep	? 🐒
EasyLift	Beam On - 4	n Current: • 50 pA	-
Direct adjustments	High Voltage 📃 🕇	2.00 kV	•
Sample preparation	Beam		? 🐒
Column presets	Stigmator	Beam Shift	
Alignements	Scan Rotation		? 🐇

Experiment

Click in the top left quadrant (the SEM window). Then click the top icon (Beam control) - Click Wake up

- Click beam on

Under the column tab, click beam on. You should hear a click from the back.

Stage navigation



Stage		?	Goto / stop buttop stage
Actual	-	Go To	
🗌 X	0 mm	4	movement
Y	-0.0001 mm	4	1
□ Z +		#	Stage position in X,Y, Z, tilt
T	0.0 °	▼ [4]	angle and rotation
R	0 °	▼ 4	
Comp	oucentric Rotation	Z-Y Link	
Last Pos	ition	Add	4
	-	Update	Saved positions
		Remove	
		Remove All	
Touch	Alarm Enabled		1
	1310/ J		Stage view from Nav Cam
1 and	- 1		
	° / ·	i a	
ő	• • • •	e =	
20			
P	0.000	and Lot	
	000		
	1		
1			

Demonstration: The inside of the e-beam

Prerequisites:

Sample loaded

Electron beam on

Learn about the SEM basics



Interaction volume and the effect of the acceleration voltage (=HT, HV)



D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland



Interaction volume and the effect of elemental composition

Secondary electrons vs backscattered electrons



D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland

Prereau	isites:
---------	---------

Sample loaded

Electron beam on

Learn about the working distance

	Experiment: Setup lens & detector
In the top menu:	
1. Select the Standard case	
2. Set magnification: low (< 100)	69 × •
3. Acceleration voltage: 2-5 kV	5.00 kv •
4. Beam size: 100pA – 800pA	0.40 nA
5. Select the ETD dectector in the topleft quadrant	det ETD

These settings are generic and not perfect yet.

- Unpause the e-beam image (click the green pause button in the image or F6).
- Refresh contrast brightness (Click *Tools* > auto contrast / brightness or **F9**)

Focus an image

If this is the first image: use e.g. a sample in the central slot² or the multisample holder to find an edge to focus.

Experiment: Focus your image

You can focus the image using:

- The coarse / fine buttons on the right of the panel
- OR Click + hold RMB in the ebeam image and move the mouse right and left

Important: after focussing, click the **LINK button** in the top row. $\boxed{\mathbb{A}} \rightarrow \boxed{\mathbb{A}}$ This allows the system to estimate the distance between the pole piece and the stage. You will end up (after homing) with a working distance around 60 mm (Z value in Navigation \Rightarrow stage).

² This is the reason why you always want to have at least one stub in the central slot. D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland

Demonstration: Eucentric height

Prerequisites:

Focused imaged Magnification: 20-5000X

Learn about the Eucentric height

This is one of the two key alignments. Every sample on every session on the FIB-SEM must be aligned to eucentric height. Failing to do so will damage the instrument.

The magic number is 7 mm

At the eucentric height position, the stage tilt and the beam axes intersect.



When the stage is tilted or rotated in any direction, this point remains focused and (almost) does not shift. At the eucentric position, one can use various system components in a safe and optimal way (e.g. GIS, Ion beam, EasyLift).

Eucentric position should be adjusted after loading any new sample

Experiment: Eucentric height position

Presettings

- 1. Assuming you homed the stage during pumping
- 2. Assuming you focused the sample or stage, and the stage is linked 🖉 (see p. 18)
- 3. Get a live Chamber View (bottom right): 5KV, Mag about 500-1000X, 0.4 nA
- 4. Click the navigation button on the right side of the screen 🚸

Procedure: decrease $Z \rightarrow adjust$ focus $\rightarrow link$ stage

- Usually, the initial Z will be somewhere 60-90 mm in a homed stage. Go in steps of 20-30 mm towards 20 mm
- E.g. Starting Z = 60mm \rightarrow Focus image, click link stage.

Set Z to 30 mm in 🚸 → Focus, link,

Set Z to 20 mm \rightarrow Focus, link,

Set Z to 10 mm \rightarrow Focus, link, 7 mm

Eucentric heigth (7 mm) reached!

- The exact starting height is dependent on the height of your sample. The values here are for a flat $\frac{1}{2}$ inch stub.

- Be conservative! Assure you are never going to touch the polepiece!!!

- When moving, Keep the mouse pointer on the stop button (-> stage) and make sure the chamber view is live.

- Clicking the Link stage button updates the working distance.
 - When the stage height is set to 7.00 mm:
 Stage > Z=7.00 mm
 - Set the eucentric height: press CTRL + f (Acknowledges a working distance of 7 mm)
 - Optional: Double click LMB in the chamber view to update the 7 mm marker.

Demonstration: Nav Cam photo

Prerequisites:

Focused imaged Eucentric height set

Learn about the the Navigation camera

The navigation camera (Nav Cam) will show the stage from a bird's eye perspective and can be used as an interactive map. It is linked to the positions and can be moved by double-clicking. CTRL + / - will digitally zoom in / out.

Moving around in a digitally zoomed nav cam picture (see p 24):



Experiment: Take Nav Cam Photo

A Nav Cam photo must be taken after proper eucentric height setup³ because the NavCam has a focal point that is reached when the stage is linked and in eucentric heigth.

Stage > take Nav Cam Photo

The green cross marks your current position.

The yellow cross (if available) marks the centre of the quadrant (View > Centre cross, or SHIFT + F5).

³ Which is why it is not recommended to do this during the pumping of the chamber (see p. 12) D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland

Prerequisites:

Eucentric height

Nav cam recorded

Lens alignment, Beam shift reset, astigmatism, Videoscope, ...

Stage module



Three modes are possible via the list box:

• Actual mode (default) – shows actual position coordinates in the edit boxes.

• Target mode – activates when clicking on a stored position or when editing a coordinate value.

• Relative mode – used to move stage by a given value and to repeat it several times if needed.

Clicking on the Go To button drives the stage to a new location. During the stage motion the Go To button changes to the **Stop button, which stops the stage immediately.**

If the symbol next to Z is ⁺⁺: You must first **LINK THE STAGE**!

Demonstration: Basic SEM imaging

Experiment: Moving around in X and Y

LMB double click in the e-beam image will move the stage. LMB double click in the Nav Cam image will move the stage. \rightarrow , \leftarrow , ... will move the stage roughly ½ field of view.

Location list

The Location list shows stored stage positions. There is the Last Position (the stage position before latest movement) in the list by default, which is updated during stage operations. The position selected becomes the actual active position and it is highlighted in the list and also in the map area. Clicking on a position name allows a user to edit it.

Map area

In the map area, the stage schema is represented showing all stored positions, which are listed in the Location list.

Reduced area (=Focus window)

This mode is useful when focusing and correcting astigmatism as the imaging update is faster in the smaller area.



Experiment: Reduced area (F7)

What is it? Reduced area allows to set and select a sub-area in the image. This assures a faster refreshing rate and is useful to focus, stigmate, ...

In case of electron sensitive samples, alignments can be made in the reduced area, away from the region of interest.

Beam shift reset



Controls the beam shift with respect to the objective lens axis. It is useful for fine imaging shifts without stage movement.

Right-clicking on the Beam Shift 2D box opens a contex	t
menu:	

• Zero sets the Beam Shift value to zero

• **Reset** sets the Beam Shift value to zero and moves the stage to compensate the resulting imaging shift (same as the Stage menu / Beam Shift Reset function).

	Coarse
1	Normal
	Fine
	Adaptive Sensitivity
	Zero
	Reset

Experiment: Reset beam shift

What is it? Beam shift is often needed for fine alignments: it allows the image to be moved very precisely (much more precisely than a piezo stage would do). Resetting the beam shift at the start of your session is recommended to have the full range of the beam shift available.

Beam control > Beam > RBM click > Zero

Digital Zoom

RAM based navigation across enlarged views. Can be used on all quadrants and detectors.

Experiment: Digital zoom

Click on the + / - button to enlarge / reduce the view in a selected display or select a zoom factor from the drop-down list.

You can use CTRL + or CTRL - on any quadrant to digitally zoom in. To position the digital zoom, select the detectors menu on the right side of the screen and in the digital zoom tab, grab and hold the green box to move it around.

Contrast and brightness can be adjusted, too. These double for the image (Contrast and Brightness buttons on the physical USB panel)

Detectors	? 🗩
Contrast -+ 61.69 %	
Brightness - + 41.02 %	
Digital Zoom	?
- + 1.0 × ·	
	and the second sec

When the digital zoom is applied, the magnifying glass icon appears in the appropriate display.

Enhanced image module

Various digital image enhancements applied to the active display independently. In case a user changes the default settings of LUT / Color / Process tab, its background changes to orange. The digital processing can be applied to any live, paused or loaded image, including an optical one.

LUT Mix 3 Mix 4 Color Process ?	LUT Mix 3 Mix 4 Color Process ?	LUT Mix 3 Mix 4 Color Process ?
D. Contrast 1.0 D. Brightness 0.00	1+2 1+2+3	•
	Gray 16 bit	
Gamma 1.0	Source 1 Source 2	
	Invert Invert	
	Source 3	
	Invert Save	Save

D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland

Lens alignment



Lens alignment

- Lens alignment is needed at magnifications > 2000X. It assures that there is no beam shift during focusing.

- Select a reduced area, preferable with sufficient contrast (press F9)

- Exposure time < 500 ns
- Column: standard, 2-5kV, 0.40 nA.

- From the top menu: Start the lens alignment. The goal is minimal movement of the object in the reduced area

- Grab the green cross that appears in the center



- Move the cross with the LMB until the movement of the object is minimal
- A slight rotation is not problematic. Unclick Lens alignment, reduced area

Astigmatism

Astigmatism in an electron magnetic lens refers to an optical aberration where the lens does not focus the electron beam to a single point along the optic axis. Instead, electrons passing through different meridional planes (typically perpendicular to each other) are brought to different focal points along the optical axis. This occurs because the lens exhibits different focal lengths for rays in orthogonal directions, often due to imperfections or asymmetries in the magnetic field or the physical construction of the lens.



Astigmatism

To correct for astigmatism (typically at magnifications > 10 000X):

- Find an object, preferably round
- Make a reduced area
- Use the X and Y under stigmator on the panel OR Shift + RMB and hold. Move left right
- / up down to correct
- AutoStigmator. First make a reduced area, then use the auto stigmator
- Switch to FFT to manually correct the astigmatism.

Videoscope

This tool facilitates manual contrast and brightness optimisation. Three yellow horizontal lines are placed over the display:

Videoscope

- Topline: whitest pixel
- Middle line: median grey
- Bottom line: blackest pixel

The oscillogram signal amplitude reflects the contrast / brightness of the just scanned line. If the oscillogram is cut by the bottom / top line, the signal level is clipped in black / white, Clipping should be avoided.

1. Select a slow scan in an active display.

2. Activate the Videoscope in an active display (F3 / clicking on toolbar icon / Scan menu) or in all live displays (Shift + clicking on toolbar icon).

3. Reduce the contrast to zero (front panel) and adjust the brightness level to the lower dashed line (black).

4. Increase the contrast so that the signal level just touches the upper dashed line (white).

5. If necessary, adjust the brightness level so that the average signal level is roughly in the middle.

6. The high and low peaks should just touch the dashed lines.

Note: Auto Contrast Brightness (F9) usually corrects these settings automatically.



Sets the magnification for the SEM image in the upper left quadrant.

High voltage (=acceleration voltage, =high tension)

Defines the interaction volume: Higher KV = less surface information, better resolution



Is the flux of electrons in the beam (electrons per second per area).

Higher beam current = better signal, more beam damage, lower resolution. Unlike the Tescan SEM, the beam current at the FIB SEM is calibrated (units in pA to nA).

- + 1 μs

Demonstration: Imag	ing setting	S
----------------------------	-------------	---

Prerequisites:

Eucentric height

Nav cam recorded

Save optimization and image settings

Image setup and filters

The image scan is a combination of:

Image size (in pixels). The larger, the longer scan will take.

Dwell time the amount of time each pixel is scanned.

Scan interlacing Splits an imaging area into n number of blocks. The first line of each block is scanned followed by the second one etc. This imaging method significantly **reduces sample charging**. Abbreviation: SI

Line integration Each line scan is repeated n times. Data are integrated.

This imaging method reduces sample charging

and improves overall image quality. Abbreviation: LI

The use of these tools is visible in the dwell time adjuster:

Image averaging

Three settings are possible:



Live. The default mode. One frame follows the other.



Average. Continuously averages n images to improve the signal to noise. Setting n > 4 will create difficulties to operate a live image (focusing, moving,

...).



Integrate. Collects n images and integrates (=sums) them. Not compatible with live imaging.

D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland



• Snapshots

A snapshot (**CTRL+F4**) is a quick recording using preset choices. The preset can be changed in the top menu or changed ad hoc using the dropdown menu. The settings for the shortcut can be set in the preferences:

Click the top left quadrant (SEM) \rightarrow Menu Tools > preferences > Scanning

\$	Electron Beam					12	1024	2048	4096	42	884	1768	3536
	25 ns	Snapshot Preset (F4)		•		88 × 5	36 ×	772 ×	× 44	2×4	24 ×)48 ×	× 96(
	50 ns	Settings				76	Ħ	200	61	2	10	20	4
\odot	100 ns	Dwell Time	300 ns		50 ns								
	200 ns	Line Integration	1		100 ns								
	300 ns	Resolution	768 × 512	-	200 ps								
	500	Line Time	284.10 μs		200 115								
	500 ms	Frame Time	147 ms		300 ns								
	1 µs	Refresh Rate	6.82 Hz		500 pr	:::::				:::			
	3 µs	Integrate	1	-	500 115								
9	5 µs	Bit Depth	16 bit	-	1 µs					:::			
	10 us	Drift Correction			3 us								
	20	Continuous Scan											
	20 µs	Action	None	-	5 µs								
	30 µs	Display	None		10 us					:::			
0	300 ns		Savo										
0	10 µs		Save As		20 µs								
			Save AS	Derauit	30 µs	:::::				:::			

The presets allow to set 8 bit (**don't!**) / 16 bit (**do!**) settings, and the proceeding function, e.g. Save as (with a dialog for a file name) or Save (with a generic name).

```
Photo
```

A photo is a high quality, slow recording using preset values. The preset can be setup as explained above for the acquisition of a photo (**F2**). There is no *ad hoc* change of dwell times / image sizes.

Demonstration: Image presets

Prerequisites:

Eucentric height

Nav cam recorded

Image parameters sets

Scanning presets simplify switching imaging conditions

By default, there are 6 factory toolbar Scanning presets (labelled s#) for each beam. Clicking on any one starts image acquiring with corresponding parameters.

Scanning presets (s1 – s6) s1 s2 s3 s4 s5 s6

At the top right, 6 scanning presets can be chosen and edited. Actions:

Select a preset: LMB

Edit a preset: RMB > edit... (click apply to save)

	Activate	Click	Scanning presets		×			
	Multiple Activate	Shift + Click	Basic Setup					
	Multiple Activate	SHILL + CILK	Name	LiSlowScan_00	1			
	Snapshot	Ctrl + Click	Resolution	768 × 512	•			
	Multiple Conservation	Ctrl + Shift + Click	Dwell Time	300 ns	•			
	Multiple snapshot	Ctri + Shirt + Cilck	Bit Depth	8 bit	-			
☆	Update with Current Settings		Filter Setup					
			Scan Interlace	1				
	Edit		Line Integration	1				
	Import Scopping Prosets		Frame Average	8	•			
	import scanning Presets		Frame Time					
	Export Scanning Presets		Image Acquisition					
	Reset to Default		Integrate	1	•			
	Reset to Belduit		Drift Correction	No	•			
			Continuous Scan	No	•			
			Action	Save As	•			
			Acquisition Time	968 ms				
			Shared Settings					
			Mains Lock	No	~			
			Start scan on left click	Yes	•			
				CI	ose			

Mains Lock When ticked, the scanning (line sawtooth signal) is synchronized with the mains AC oscillation. This greatly diminishes blurring and jittering of the electron imaging resulting in smooth image edges at higher magnifications.

D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland

Demonstration: Align Feature / scan rotation

O Horizontal

OK

Vertical

Prerequisites:

Eucentric height

Nav cam recorded

Image parameters sets

Rotate the scan angle and align your sample of interest

The scan rotation can be adjusted to align the image with possible objects.

Experiment: scan rotation Click the Align feature button in the top row. A window will appear: - You can now by LMB click and draw a yellow dotted line in the electron beam quadrant. The angle of the line will be shown. You can now choose to: Rotate the stage Align Feature Rotate the scan Move points or line by dragging them Position 1 Position 2 X = -26.281 mm X = -26.019 mm Y = -13.772 mm Y = -13.785 mm Click OK: the image will rotate (either Orientation Angle by scan or stage) to align the feature

3 °

Demonstration: Column cases

Prerequisites:

Eucentric height Nav cam recorded

Image parameters sets

Learn about the three column cases: Standard, OptiPlan, OptiTilt

The electron column can be operated in different Use Cases optimized for specific applications: Standard, OptiTilt and OptiPlan.



In the OptiTilt and OptiPlan Use Cases: primary electrons are accelerated by the potential of the Acceleration tube (A-Tube) and pass through the column at high energy (i.e. reducing aberrations). They are decelerated between the T1 detector and Secondary the sample. as well as backscattered electrons are also collimated the final by the A-Tube into lens electrostatic field and detected by detectors T1 and T2.

• Standard mode

This is the basic mode, used for survey and TEM sample preparation mode. It is ideal for navigating and reviewing sites at lower magnifications. Maximum e-beam current is 13 nA.

- A-tube: deactivated
- Typical detector: ETD
- Advantage: largest field of view, all currents available
- Disadvantage: not suitable for high resolution (no beam deceleration)

* OptiPlan mode

This mode is used for ultrahigh resolution electron imaging at 0° of the sample at short working distances (1 - 2 mm). Exceptionally not at 7 mm). The A-tube is at the highest potential and the T1 and T2 detectors should be used. Full range of the Beam Deceleration mode (see below) is available.

- Typical detector: T1 & T2 (note that the ETD detector signal will drop)
- WD can be set to < 5 mm (you anyway will not tilt, it is only useful at 0°)
- Advantage: high resolution, more signal
- Disadvantage: smaller field of view, not all currents available

* OptiTilt mode

The OptiTilt Use Case is optimized for operation at eucentric working distance (i.e. 7 mm, 52° tilt). It is ideal for FIB cross section preparation and subsequent ultra-high resolution electron imaging at low accelerating voltages. The A-Tube is on the high potential, the T2 or ETD detectors are used to obtain topographical images and T1 to gain the

composite information. A small positive stage bias (see below) to improve the image quality is applied automatically when the stage is tilted.

- Typical detector: T1 & T2 (note that the ETD detector signal will drop)
- WD should be at 7 mm (or about there, depending on BCP)
- Advantage: high resolution at 52°
- Disadvantage: smaller field of view, not all currents available

Demonstration: Detectors

Prerequisites:

Imaging of samples

Overview of all installed detectors

The system has 8 detectors installed.

*	-
4	
۲	
ww	
3	
tlt	

STANDARD DETECTOR: ETD (=Everhart Thornley detector = standard SE detector)

Detects low energy (=secondary electrons) electrons generated by the primary beam interaction with the specimen surface. It is permanently mounted in the chamber

over and to one side of the sample.

De	tector ETD
	-
Mode	Secondary Electrons
Grid voltage	- + 250 V

Mode: keep this the Secondary electrons **Grid voltage**: Bias applied by the collector. The higher, the stronger SE (and noise) are attracted. Sweet spot at +250 V. Negative voltage repels SE from the ETD detector and only BSE are detected.

IN LENS BSE DETECTOR: Trinity 1

Electrons generated by a primary beam can be collected by the in-lens Trinity

detectors T1 and T2, which are located inside the final lens. Whenever the T1 or T2 is selected, the CCD camera infra-red LED's are switched off not to emit the photons.

The T1 detector is primarily designed to **collect backscattered electrons** (BSE) and provides composite sample contrast. In the OptiTilt and

Detector Settings				
	Detector T1	-		
Segment	t A			
Segment	t B			
A + B				
O A - B				

OptiPlan column Use cases, it detects backscattered electrons through the whole range of accelerating voltages.

In the Standard column Use case, the T1 provides strong BSE contrast at accelerating voltages of 5 kV and higher.

The active detector area is split into two halves and the detector can be operated in four modes. Apart from the composite mode A + B the detector can be operated in the topographical mode A - B, where a pseudo-topographical imaging with suppressed atomic number contrast and maximum topographical contrast is obtained. Signal from each half can be also collected separately in mode Segment A / B.

IN LENS SE DETECTOR: Trinity 2

The T2 detector is primarily designed to collect secondary electrons (SE) and provides information of the sample topography in the OptiTilt and OptiPlan column Use cases.

It can be operated in Standard column Use case as well. In this case, backscattered electrons are collected, but the accelerating voltage must be 5 kV at least. The intensity



of the signal in the Standard column Use case is dependent on the beam current and working distance – the higher the beam current and shorter the WD the higher the T2 signal. There are no options to select.

RETRACTABLE BSE DETECTORS: CBS and ABS⁴ detector

The CBS uses concentric segmentation of the detector diode to distinguish between



backscattered electrons scattered close to the beam axis – inner segment (preferentially composite contrast) and electrons scattered far from the beam axis – outer segment (more topographical signal).

⁴ Concentric and Annular backscattered electron detector

D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland

Do not use the detector with high ion beam currents (the solid-state diode can degrade). When working with ion beam for longer time, retract the detector! Be aware of sample and final lens collision when tilting large samples!

Select the CBS or ABS from the Detector Settings module > Detector list box. Choose the required diode segment(s) by selecting the relevant radio button. The Custom mode is used to define the segments to be used for detecting. Clicking on the + / - sign in particular segment activates it to add (yellow background) / subtract (blue background) the segment signal. By double-clicking on the segment its background turns grey and it is switched off. The Contrast button equalizes signals (contrast) from different segments not to override one another. Distribution of electrons collected by detector segments changes with setting of working distance, lens mode and Beam Deceleration mode. It is also possible to set different concentric segments in particular displays and thereafter to use the Enhanced Image module / Mix 3 or Mix 4 tab to mix color coded signals to create color images.

SCANNING TRANSMISSION ELECTRON MICROSCOPY: STEM and STEM+ detector

This detector is retractable and is mounted below the sample. It is only useable in conjunction with the rowbars. The detector has selectable segments that allow to operate the diode in

Bright Field, Dark field and High Angle Annular Dark Field mode. The resolution of the STEM is 0.9 nm maximum.





EDX DETECTOR: analytics

Not part of this introduction.

Demonstration: Direct adjustments – Source tilt

Prerequisites:

Poor signal conditions in the e-beam

Fine-tuning of the e-beam geometry

Notice: only perform the here-mentioned procedure, and only when it appears to be necessary. Do not perform any other direct alignments or click any other buttons (see <u>p 4</u>).

WHEN: if all other alignments (eucentric height, camera settings, beam current, ...) appear correct, but the e-beam is providing a very noisy, low signal.

This algorithm does not affect the ion beam!



A green X marks the centre. When finished, the be concentric around the green X.

Press F9 to correct brightness contrast.



Experiment: Source tilt alignment

D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland

Demonstration: Beam deceleration

Prerequisites:

Imaging of samples

Learn about beam deceleration

The Beam Deceleration (BD) mode is based on a negative voltage (bias) applied to a stage (i.e. a sample). The electrical field between the sample and the nearest surface over (a column bottom or a detector) is formed, acting as the additional electrostatic lens. Its power is described by the Immersion Ratio parameter:



As the sample is at the negative potential with relation to the ground and detectors, the initial SE and BSE energy (when leaving the surface) is accelerated by the Stage Bias before the detection.

The higher the Immersion Ratio, the lower is the difference between SE and BSE energies when detected. Signal electrons are accelerated upwards and deflected towards the column axis.

The secondary electrons

The SE's have a low initial speed and they are usually absorbed into the detector central hole, continue through and can be detected by in-column detectors (T2). Equally like the BSE heading upright.

The back scattered electrons

Conversely the BSE heading nearly parallel to a surface (which normally cannot be detected) are driven to a detector.

By changing the Stage Bias an output angle of electrons leaving a surface can be changed.

OptiPlan and OptiTilt mode will apply a stage bias.

The stage bias, typically 20V, will create the immersion effect. This increases the electron yield on the in-beam detectors T1 and T2. When active, note the small arrow after the KV value. The beam deceleration accelerates the beam in the column by 8 kV (this reduces chromatic aberration in the lenses) and decelerated before the beam exits the pole piece.

Beam Deceleratio	n		?		
On	Stage	Bias -50.00 V	· ·	HV N 6.00 KV	bias 3.00 kV



Experiment: Shutdown

1. Electron beam and ion beam: same procedure as standby (see above)

2. Sleep

Click the sleep button in the System menu, just above the Column window.

System	?	System	?
Wake Up	Sleep	Wake Up	Sleep
Column	? 🏀 Beam Current	Column Beam C	? 🍪
Beam On	-+ 0.50 nA	Beam On -+ 0	.50 nA
High Voltage	4 II → /	High Voltage 🖣	

The green bar below the Beam On button of the ion beam turns yellow, then red and reduces in size.

Summary Login • Vent the chamber, then load your samples (tallest on the right) • Pump the chamber + Assure the door is tightly closed / Lean against the chamber door Start •Home the stage during pumping Switch on the beams Choose a proper HV setting / Beam intensity • go to minimum magnification (< 100X), find your sample Beam Focus and link your image • Eucentric height: decrease the Z position iteratively until Z=7mm. CTRL+F • Perform beam coincidence point: go in steps to 52°. Use MMB to shift the Z position • Match e-beam and ion beam using Shift + LMB. Use ion beam shift. Sample Define image scanning properties Record your data Save data on D:/sharedData Data Secure your data through the support PC • Remove your samples and pump the chamber + + Assure the door is tightly closed / Lean against the chamber door End Switch off the e beam and the ion beam and Log out

D Vanhecke | Adolphe Merkle Institute | University of Fribourg | Switzerland

adolphe merkle institute UNIVERSITY excellence in pure and applied nanoscience SWITZERLAND